


Alginate Methacryloyl (AlgMA)

Product component

Item	Character	Package Size	Notes
A: AlgMA	White spongy	0.2 or 0.5g/bottle	Keep in dark
B: Photoinitiator LAP	White powder	0.025g/bottle	

This instruction applies to EFL-AlgMA-50K/300K

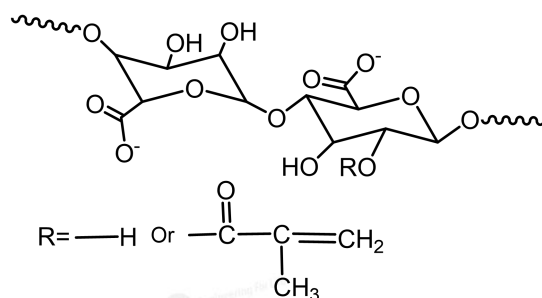

AlgMA molecular structure

product introduction

Alginate methacryloyl (AlgMA) is a double-bonded modified sodium alginate, which can be cross-linked and cured into the gel through UV and visible light in the presence of the photoinitiator. Compared with traditional divalent ions (calcium ions, etc.) crosslinking, AlgMA photo-crosslinked method is highly portable and has good internal homogeneity in the gel. AlgMA light-curing hydrogels have a three-dimensional (3D) structure suitable for cell growth and differentiation, and both -OH and -COOH in the structural units can be used as active sites for chemical reactions. In addition, AlgMA hydrogel also has good mechanical properties. And the 3D hydrogel micro-scaffolds constructed by it have adjustable mechanical and chemical properties.

Applications

3D cell culture, biological 3D printing, tissue engineering, etc.

Storage

Dry kit: room temperature, 3 months; 4°C, 12 months; -20°C, 18 months. **Sterile solution:** 4°C (in dark), 7 days; -20°C (in dark), 6 months. **Please note that repeated freezing and thawing of the solution will affect the performance of the product, so it is best to prepare it when using it.**

Period of validity



企业微信公众号
扫描右侧二维码
获取更多信息

The date of manufacture is shown in the package.

Solution preparation

1. Prepare 0.25% (w/v) standard solution of initiator

- (1) Add 10ml PBS into the brown bottle containing initiator LAP (containing 0.025g LAP);
- (2) Heat and dissolve the solution in a water bath at 40-50°C for 15 minutes, shaking several times.

The LAP standard solution can be stored for 12 months at 4°C in dark.

2. Prepare AlgMA solution (5-10% (w/v) is recommended for AlgMA-50K, 1-2% (w/v) is recommended for AlgMA-300K)

- (1) Take the required mass of AlgMA into the centrifugal tube;
- (2) Take the required volume of initiator standard solution and add it to the above container;
- (3) Stir and dissolve at room temperature in the dark for 0.5-1h;
 - The viscosity of the AlgMA-300K solution is large. The dissolution time can be extended appropriately. Pay attention to seal to prevent moisture volatilization;
 - It is recommended to use centrifugal machines to remove bubbles from the system (3000~5000rpm, 2-3min);
- (4) Sterilize the AlgMA solution with a 0.22μm sterile needle filter and keep in dark.

Suggestions for 2D cell culture

- Inject AlgMA solution into the well plate;
(96-well plate: 50-100 μL / well, 48-well plate: 100-300 μL / well, 24-well plate: 300-500 μL / well);
- Irradiate the wells with 405nm light for 10-30 seconds to gelate, the gel strength can be adjusted by the time and intensity of the light;
- Add medium to the wells. Place the plate in a 37°C incubator for 5 minutes. And then wash the sample and remove the medium;
- Add the cell suspension to the well plate. Change medium, observe, and photograph according to experimental design. (No special requirements for operation procedures).



企业微信公众号
扫描右侧二维码
获取更多信息

Suggestions for 3D cell culture

- Cells are collected and resuspended in pre-warmed AlgMA solution to prepare the cell suspension;
- Add cell suspension into the well plates;
(96-well plate: 50-100 μ L/ well, 48-well plate: 100-300 μ L/ well, 24-well plate: 300-500 μ L/ well)
- Irradiate the wells with 405nm light for 10-30 seconds to gelate, the gel strength can be adjusted by the time and intensity of the light;
- Add medium to the wells. Place the plate in a 37°C incubator for 5 minutes. And then wash the sample and remove the medium;
- Add fresh medium and incubate for a long time. Change medium, observe, and photograph according to experimental design. (No special requirements for operation procedures).

Tips: Do not look directly at the light source.



企业微信公众号
扫描右侧二维码
获取更多信息